



METALS SAMPLING INSTRUCTIONS (IEPA)

Metals - contain the following:

- ▶ **1 – 250ml plastic bottle preserved with Nitric Acid**
“DO NOT EMPTY OR RINSE OUT PRESERVATIVE”

- 1) Please fill bottle to the neck from the same location listed on the COC.**
- 2) Identify sampler name, date and time of collection, and relinquished by on the COC.**
- 3) If you get any acid on your skin, flush with water for several minutes. If necessary call your physician.**
- 4) Metals do not have a short hold time and do NOT require expedited shipping. No regulatory cooling or refrigeration required.**

NO ICE IS REQUIRED!!!!